



FH5
Forced Air Oven
Operation and Service Manual

TestEquity Inc.
2450 Turquoise Circle
Thousand Oaks, CA 91320

Support: 877-512-3457 Toll Free
805-480-3697

Corporate: 800-732-3457
805-498-9933

<http://www.testequity.com>

Table of Contents

Chapter 1 – Safety Instructions	1-1
Introduction	1-1
Safety Notices	1-1
Chapter 2 – Receiving and Installation	2-1
Inspection	2-1
Return Shipment	2-1
Installation	2-1
Power Requirements	2-1
Power Wiring	2-2
Location	2-2
Lifting / Handling	2-2
Leveling	2-2
Initial Burn-In	2-2
Initial Cleaning	2-2
Chapter 3 – Operation	3-1
Front Panel Controls	3-1
Power Switch	3-1
Overtemperature Protection Thermostat	3-1
Overtemperature Light	3-1
Temperature Controller	3-1
Intake and Exhaust Knobs	3-1
Operating the Oven	3-2
Turning the Oven On	3-2
Setting the Temperature Controller	3-2
Setting the Overtemp Protection Thermostat	3-2
Convenience Outlet	3-2
Chapter 4 – Temperature Controller	4-1
Introduction	4-1
Single Set Point Operation	4-1
Single Set Point Operation	4-2
Programmed Operation	4-2
Chapter 5 – Maintenance	5-1
Cleaning	5-1
Storage	5-1
Calibration	5-1
Troubleshooting Guide	5-2
Parts List	5-5
Specifications	5-5
Wiring Diagram	5-6
Chapter 6 - Warranty	6-1

Chapter 1 – Safety Instructions

Introduction

Follow all CAUTION notices to prevent damage to the chamber or your test sample. Failure to follow all CAUTION notices may void your warranty. CAUTION may also indicate a potentially hazardous situation which, if not avoided, may result in minor or moderate personal injury.

WARNING indicates a potentially hazardous situation which, if not avoided, could result in death or serious injury.

The safety alert symbol  precedes a general CAUTION or WARNING statement.

The electrical hazard symbol  precedes an electric shock hazard CAUTION or WARNING statement.

Safety Notices

 **WARNING:** These units are general purpose air ovens for professional, industrial or educational use where the preparation or testing of materials is done at approximately atmospheric pressure and no flammable, volatile or combustible materials are being heated. These units are not intended for hazardous or household locations or use.

 **WARNING:** THIS OVEN IS NOT DESIGNED TO HANDLE COMBUSTIBLE GASSES AND IS NOT AN EXPLOSION PROOF UNIT. Do not place explosive, combustible, or flammable materials into the chamber.

 **WARNING:** Some outgassed byproducts may be hazardous or unpleasant to operating personnel. If this is the case, the oven exhaust should be positively ventilated to the outside according to local regulations. A power exhaust is optionally available which greatly helps under these applications.

 **WARNING:** Do not operate near noxious fumes.

 **CAUTION:** Do not place sealed or filled containers in the oven chamber.

 **CAUTION:** If a mercury thermometer is used for calibration and breakage should occur, all spilled mercury MUST be completely removed from the chamber before continuing operation.

 **CAUTION:** Be sure that the power supply is of the same voltage as specified.

 **WARNING:** Do not cut or remove the ground prong from the power cord (if one is used). Do not use a 2-prong adapter plug.

Chapter 1 – Safety Instructions

- ⚠ WARNING:** Disconnect the unit from its electrical source before proceeding to make any electrical repairs or replacements.
- ⚠ CAUTION:** This oven is NOT designed for the use in Class I, II, or III locations as defined by the National Electric Code.
- ⚠ WARNING:** This oven is not intended, nor can it be used, as a patient connected device.

Chapter 2 – Receiving and Installation

Your satisfaction and safety require a complete understanding of this unit. Read the instructions thoroughly and be sure that all users are given adequate training before attempting to use this unit. Note: This equipment must be used only for its intended purpose; any alterations or modifications will void your warranty.

Inspection

The carrier, when accepting shipment, also accepts responsibility for safe delivery and is liable for loss or damage claims. On delivery inspect for visible exterior damage, note and describe on the freight bill any damage found and enter your claim on the form supplied by the carrier.

Inspect for concealed loss or damage on the unit itself, both interior and exterior. If any, the carrier will arrange for official inspection to substantiate your claim.

Verify that your accessory package is complete. All units should have a set of four (4) Leveling feet, two (2) shelves and four (4) shelf slides.

Return Shipment

Save the shipping crate and all packaging material until you are sure the oven has not been damaged and is functioning correctly. If for any reason you must return the unit, contact your customer support representative for authorization and supply data plate information. Make sure to include the model and serial number. The service representatives will furnish you with a return authorization number and address for return. Note: Make sure this return authorization number appears on the unit packaging and shipping papers. Units returned without proper authorization may not be accepted at the factory. For information on where to contact customer support please see the manual cover.

Installation

Local city, county or other ordinances may govern the use of this equipment. If you have any questions about local requirements, please contact the appropriate local agency. Installation may be performed by the end user.

This unit is intended for use indoors, at room temperatures between 5° to 40°C (41° to 104°F), at no greater than 80% Relative Humidity (at 25°C/77°F) and with a supply voltage that does not vary by more than $\pm 10\%$. Customer support should be contacted for operating conditions outside these limits.

Power Requirements

The power requirements are listed on the oven's data plate located on the rear. Make sure your power supply matches that shown on the data plate. **VOLTAGE SHOULD NOT VARY MORE THAN $\pm 10\%$ FROM THE DATAPLATE RATING.** These units are intended for 50/60 Hz application. A separate circuit is recommended to preclude loss of product due to overloading or circuit failure.

Power Wiring

Installation requires hard wiring and should be performed by a qualified electrical technician. The next higher circuit breaker value above the data plate amperage may be used, provided the requirements in article 422 of the National Electric Code are met or, the applicable code that applies in the country where this unit will be operating.

Location

Select a site for the oven which is free from extreme heat, cold or excessive air movement such as areas near steam radiators, stoves, other ovens, autoclaves, direct sun, heating and cooling ducts, etc. Avoid high traffic areas which may reduce the accessibility to the oven. Allow at least 2 inches of space between the oven and surrounding walls or partitions which might prevent free airflow.

Lifting / Handling

These ovens are heavy and care should be taken to use appropriate lifting devices that are sufficiently rated for these loads. Ovens should only be lifted from their bottom surfaces. Doors, handles and knobs are not adequate lifting or stabilization. The oven should be completely restrained from tipping during lifting or transport. All moving parts, such as shelves and trays should be removed and doors need to be positively locked in the closed position during transfer to prevent shifting and damage.

Leveling

The oven must sit level and solidly. Leveling feet (supplied) are to be installed in the holes at the base of the oven. Turn them counterclockwise to raise the level and clockwise to lower the level. If the oven must be transported, turn the leveling feet in all the way to prevent damage.

Initial Burn-In

It is recommended that the unit go through a “burn-in” process prior to operation. This is to eliminate the smoking of protective coatings on the element. Read the operating instructions carefully to understand the operating requirements. To burn-in, turn the OVERTEMP PROTECTION to maximum and set the TEMPERATURE controller to 200°C. Open the exhaust vent damper on the top of the oven. Run for a minimum of one hour under ventilation until smoke dissipates.

Initial Cleaning

The oven's interior was cleaned at the factory. If you need to clean it further, use a cleaner that is suitable for your application. Make sure to rinse the cleaned surface with a damp cloth, using water only, and dry the surfaces with a clean cloth. DO NOT USE chlorine-based bleaches or abrasives as they will damage the stainless steel surface. DO NOT USE spray cleaners that might leak through openings and cracks and get on electrical parts or that may contain solvents that will harm the coatings. A similar periodic cleaning is recommended.

⚠ WARNING: Never clean the unit with alcohol or flammable cleaners with the unit connected to the electrical supply. Always disconnect the unit from the electrical service when cleaning and assure all volatile or flammable cleaners are evaporated and dry before reattaching the unit to the power supply.

Chapter 3 – Operation

Front Panel Controls

Power Switch

The POWER switch controls all power to the oven. It must be in the I (ON) position before any systems are operational. The green light in the switch will be lit when the switch is ON.

Overtemperature Protection Thermostat

The OVERTEMP PROTECTION thermostat is an independent controller to protect against any failure which would allow temperature to rise past the TEMPERATURE controller's set point. The adjustment knob points to a relative scale from 0 to 10.

Overtemperature Light

The OVERTEMP PROTECTION pilot lamp will come on when the OVERTEMP PROTECTION is activated and has taken control of the oven. This pilot lamp should never be on under normal operating conditions.

Temperature Controller

The TEMPERATURE CONTROLLER is capable of both continuous single-setpoint and timed profile modes of operation.

Intake and Exhaust Knobs

The front-panel INTAKE and EXHUST knobs control the air vents, which are located on the top of the oven. To achieve near-ambient temperatures or to cool down from higher temperatures without opening the door, open both vents.

Operating the Oven

Assure that the electrical power source is properly configured and rated for the oven and plug the unit cord into the receptacle.

Turning the Oven On

Push the POWER switch to the I (ON) position. Push the timer switch to the O (OFF) position. The digital temperature display will indicate a temperature value. Turn the OVERTEMP PROTECTION Thermostat to its maximum clockwise position using a coin or flat head screwdriver.

Setting the Temperature Controller

When the oven is turned on the controller will show the temperature within the oven on the upper display. The lower display will show the temperature set point. To change the set point press the Up or Down keys until the desired temperature is indicated in the lower display. Details for entering and running a profile are described in Chapter 4.

Setting the Overtemp Protection Thermostat

The OVERTEMP PROTECTION should be initially set to its maximum position when stabilizing the set point temperature. Once the oven is stable at the desired set point, turn the OVERTEMP PROTECTION counterclockwise (using a coin or flat-head screwdriver) until the OVERTEMPERATURE light comes on. Next, turn the OVERTEMP PROTECTION clockwise just until the OVERTEMPERATURE light goes off. Next, turn the OVERTEMP PROTECTION two (2) minor increments on its scale past the point where the light went out. This sets the OVERTEMP PROTECTION at approximately 10°C above the current temperature set point.

Convenience Outlet

An electrical outlet is located on the rear of the oven. It provides power to an optional Power Exhaust Blower. The voltage present at this connector is the same as the oven's input voltage.

This outlet is controlled using the Event 3 function on the temperature controller.

Chapter 4 – Temperature Controller

Introduction


This chapter provides an overview on how to operate the Watlow 981 Temperature Controller. Details can be found in the “Watlow 982 Users Manual”. Note that the 981 and 982 are functionally identical.

⚠ CAUTION: The “Series 982 User’s Manual” is a general manual and is written by the manufacturer, Watlow, for a wide variety of applications and configurations. Not all features or functions are applicable.

⚠ CAUTION: The Watlow 981 Temperature Controller has been properly configured by TestEquity to match the oven’s system requirements and to perform optimally over a wide range of operating conditions. Improper modifications to these setup values can result in erratic performance and unreliable operation. Do not attempt to modify the setup values, unless you thoroughly understand what you are doing. If there is any doubt, please call TestEquity before proceeding.

The Watlow 981 Controller provides easy single set point operation and profile programming. The controller can store four profiles with up to 6-steps in each profile.

The user-interface is organized in menus. The Setup and Factory menu values are programmed at the factory and should need no adjustment under most circumstances. Do not attempt to modify these values unless you thoroughly understand what you are doing. If there is any doubt, please call TestEquity before proceeding. The Operation Menu has three sub-menus – System, PID and Program. The System and PID menus are factory preset and should not need any adjustment under normal circumstances. The Program menu is where you enter your programs.

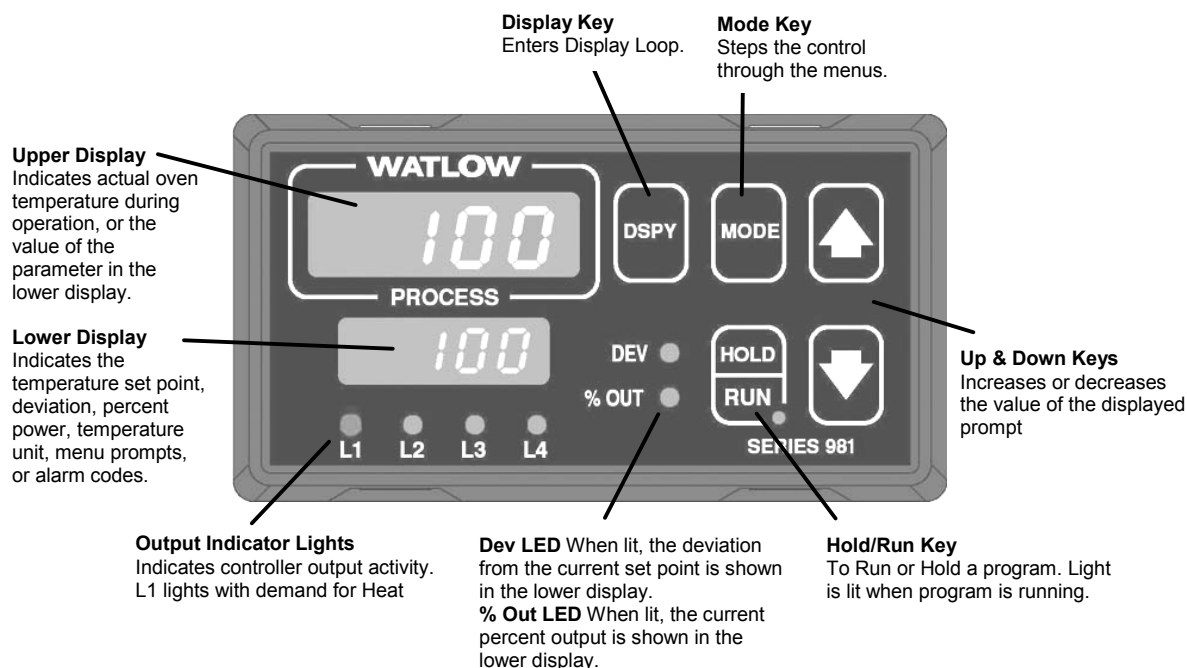

Single Set Point Operation

Normally, an average user only wants to operate the oven at one temperature. The Temperature Controller is in Single Set Point mode when it is not running a program. The Upper Display shows the actual chamber workspace temperature. The Lower Display shows the desired temperature Set Point. To enter a temperature Set Point, simply press the ▲ or ▼ key to enter the desired temperature.

Programmed Operation

The temperature controller is capable of being programmed with up to four profiles. A ramp is a controlled change in temperature over a specified period of time. A soak is a hold at particular temperature for a specified period of time with up to six steps each. Profiles can be linked for programs needing more steps up to a total of 21 ramps or soaks. Refer to Chapter 7 of the Watlow “Series 982 User’s Manual” for detailed instructions.

Sometimes a user will want to operate the oven at a given temperature for a specified period of time and then cool off. To do this, a small program must be put into the controller.

The Operation Menu has three sub-menus – System, PID and Program. The System and PID menus are factory preset and should not need any adjustment under normal circumstances. The Program menu is where you enter your programs.

The Program menu is used to input time and temperature steps into the controller. The five types of steps that can be programmed are Set Point, Soak, Jump Loop, Link File and End. A Set Point step determines the desired temperature and the time taken to get to that temperature. A Soak step determines amount of time that the temperature in the previous step is maintained. A Jump Loop step tells the program to jump to another step. A Link File step jumps to another profile. An End step ends the program.

Each programmed step consists of a series of commands. For a Set Point step commands are SP (Set Point temperature), Hours (Hours), Min (Minutes), Sec (Seconds), ENT3 (Event 3 controls the outlet for the optional power exhaust). The set point is the temperature that you want the oven to go to. The Hours, Minutes, Seconds is the time you want the oven to take to reach set point. If you want the oven to reach set point as fast as possible then set these times to zero.

A Jump Loop step has three commands – JF (Jump File), JS (Jump Start), and JC (Jump Count).

The Link File step command LFIL is used to link to another profile.

End step has only one command.

An example of a simple program would be one in which you want the oven at 90 to 100 degrees, maintain 100 degrees for 5 ½ hours and then cool down to room temperature. This would be a four step program. The first step would be Set Point step. The Set Point would be 100 degrees. The time would be programmed as zero since you want the oven to get 100 as fast as possible. The second step would be a Soak Step. The soak time would be 5 hours and 30 minutes. The third step would be a Set Point step with the set point set to zero and the time set to zero. The final step would be an End step.

Chapter 5 – Maintenance

⚠ WARNING: Disconnect the power cord from the power source before performing any service or maintenance on this unit.

No maintenance is required on the electrical components. If the unit fails to operate as specified, please see the Troubleshooting Guide, before calling for service.

Cleaning

Cleaning and decontamination are recommended on a regular basis. To prepare the unit for cleaning, remove all interior parts if assembled, such as shelves and shelf clips.

First, clean the oven interior with soap and water, rinse and let dry. To decontaminate use a solution that is appropriate for your application. **DO NOT USE** chlorine-based bleaches or abrasives as this can damage the stainless steel components. **DO NOT USE** spray cleaners that might leak through openings and cracks and get on electrical parts or that may contain solvents that will harm the coatings. Use care when cleaning the door gasket to prevent damage which could impair the positive door seal.

⚠ WARNING: Never clean the unit with alcohol or flammable cleaners with the unit connected to the electrical supply. Always disconnect the unit from the electrical service when cleaning and assure all volatile or flammable cleaners are evaporated and dry before reattaching the unit to the power supply.

Storage

To prepare the unit for storage, remove all shelves and shelf clips and disconnect the power supply. Be certain that the chamber is completely dry and door is positively secured in the closed position. See Lifting / Handling for proper transport procedures.

Calibration

It is recommended that calibration is verified with the unit is installed in its working environment, and has been stable at set point for several hours. Place a reference temperature measurement device in the oven. Be certain the sensor does not touch any shelving. Allow again for the temperature to stabilize until five (5) consecutive readings at one minute intervals show no temperature change. Compare the reading on the reference thermometer with the front panel digital display. If there is an unacceptable difference, refer to the calibration instructions in the “Series 982 User’s Manual”. If you simply want to enter an offset to compensate for uniformity in the oven (not input calibration errors), use the “Calibration Offset” function.

Troubleshooting Guide

TEMPERATURE	
Temperature too high.	1. Controller set point is too high. 2. Controller failure – call Customer Service.
Display reads error code “EX4”.	1. Probe is unplugged, is broken or wire to sensor is broken – trace wire from display to probe.
Temperature too low	1. OVERTEMP PROTECTION set too low. 2. TEMPERATURE controller set too low. 3. Unit needs time to recover from door opening – wait for display to stop changing. 4. Unit not recovered from power failure or being turned off – ovens will need several hours to warm up and stabilize. 5. Element failure – compare current draw to data plate. 6. Temperature controller failure – call Customer Service. 7. OVERTEMP PROTECTION failure – confirm with front panel lights that safety thermostat is operating correctly.
Oven will not heat over a temperature that is below set point	1. Confirm that fan is moving and that amperage and voltage match data plate – check for air movement in chamber. 2. Confirm that set point is set high enough – turn OVERTEMP PROTECTION all the way clockwise and see if OVERTEMPERATURE light comes on. 3. Check connections to sensor. 4. Check calibration using a calibrated instrument.
Unit will not heat up at all	1. Check amperage – amperage should be virtually at maximum rated (data plate) amperage. 2. Do all controller functions work? 3. Is the Safety Thermostat set high enough? For diagnostics, should be fully clockwise with the OTP light never on. 4. Has the fuse/circuit breaker blown? 5. Has timer turned unit off?
Will not maintain set point	1. Assure that set point is at least 10 degrees over ambient. 2. See if ambient is fluctuating; check for adjacent open doors or HVAC duct openings – stabilize ambient conditions.
Display and reference thermometer don’t match	1. Calibration error – Controller needs re-calibration. 2. Temperature sensor failure – call Customer Service. 3. Controller failure – call Customer Service. 4. Allow at least two hours to stabilize. 5. Verify that reference thermometer is certified.
Calibrated at one temperature, but not at another	1. This can be a normal condition when operating temperature varies widely. For maximum accuracy, calibration should be done at or as close to the set point temperature.

Chapter 5 – Maintenance

TEMPERATURE (continued)	
Indicated chamber temperature unstable	1. $\pm 0.1^{\circ}\text{C}$ variation is normal. 2. Make sure fan is working. Verify movement of air in chamber. 3. Room temperature is radically changing – either door opening or room airflow from heaters or air conditioning. Stabilize the ambient conditions. 4. This may happen if exhaust stack is 100% open or if power exhaust is cycling – adjust stack to at least $\frac{1}{4}$ closed. 5. Sensor miss-located, damaged or wires may be damaged. Check mounts for control and overtemperature sensors, then trace wires or tubing between sensors and controls. 6. Defective temperature controller. Call Customer Service. 7. OVERTEMP PROTECTION set too low. Be sure that OVERTEMP PROTECTION is set more than 5 degrees over temperature set point; check if OVERTEMPERATURE light is on continuously; turn OVERTEMP PROTECTION knob completely clockwise to see if problem solved then follow operating instructions for correct setting. 8. Electrical noise. Remove nearby sources of RFI including motors, arcing relays or radio transmitters 9. Bad connection on temperature sensor or faulty sensor. Check connectors for continuity and mechanical soundness while watching display for erratic behavior; check sensor and wiring for mechanical damage. 10. Bad connections or faulty solid state relay. Check connectors for mechanical soundness and look for corrosion around terminals or signs of arcing or other visible deterioration.

MECHANICAL	
Motor doesn't move	1. If shaft spins freely: check connections to motor and check voltage to motor. 2. If shaft rubs or is frozen, relieve binding and retest.
Motor makes noise	1. If noise is from the motor, tap the top of motor shaft with ball peen hammer. 2. If the sound gets worse, tap the other end of the shaft – avoiding touching the blower wheel. 3. If there is no change, call Customer Service.
Door not sealing	1. Adjust hinge blocks, chamber latch bracket or twist the door. 2. Confirm that unit has not been damaged and body is not square. 3. Check physical condition of gasket for tears or punctures.

Chapter 5 – Maintenance

OTHER	
Temperature is running away	1. Adjust set point to room temperature. If the unit is still heating, replace the solid state relay. 2. If cannot change any condition on the front panel, call Customer Service.
Unit or wall fuse/circuit breaker is blown	1. Check wall power source. 2. Compare current draw and compare to specs on data plate. 3. See what other loads are on the wall circuit.
Unit will not turn on	1. Check wall power source. 2. Check fuse/circuit breaker on unit or in wall. 3. See if unit is ON, e.g., fan or heater, and just controller is off. 4. Check all wiring connections, especially around the on/off switch.
Unit is smoking after initial operation	1. This is not an uncommon occurrence when first operating new units. Run at high temperature with vent open for one hour or until smoke dissipates.
Contamination in chamber	1. See cleaning procedure. 2. Develop and follow standard operating procedure for specific application; include definition of cleaning technique and maintenance schedule.
Contamination on sample	1. See “Contamination in chamber”. 2. Reduce air flow in chamber by dampening down exhaust port; be sure to verify adequate temperature uniformity at the reduced air flow. 3. Protect open samples from areas of maximum air current, e.g., inlet air ducts.


Parts List

Description	Part Number
Blower Motor	4880512
Convenience Outlet	6550515
Cooling Fan	X1000748
Cord Cap	1810502
Cord Cable	1800524
Fan Guard	2600518
Line Filter	2800503
Heating Element	9570831
High Limit Controller	1750571
High Limit Indicator Pilot Lamp	200020
Leveling Feet	200129
ON/OFF Switch	103351
Power Relay	100064
Shelf	5121154
Shelf Standards	5121189
Solid State Relay	102162
Temperature Probe	101827
Watlow 981 Temperature Controller	1750540

Specifications

Inside Volume	4.7 cu ft
Inside Dimensions*	20.5 x 20 x 20.125
Outside Dimensions*	35 x 30.25 x 38
Shipping Weight	400 lbs.
Net Weight	325 lbs.
Temperature Range	40°C above ambient to 300°C
Uniformity	±0.9°C Typical
Control Tolerance	±0.1°C Typical
Input Power	220 VAC nominal, 50/60 Hz, 12 Amps, 2200 Watts
* W x D x H (inches)	

Wiring Diagram

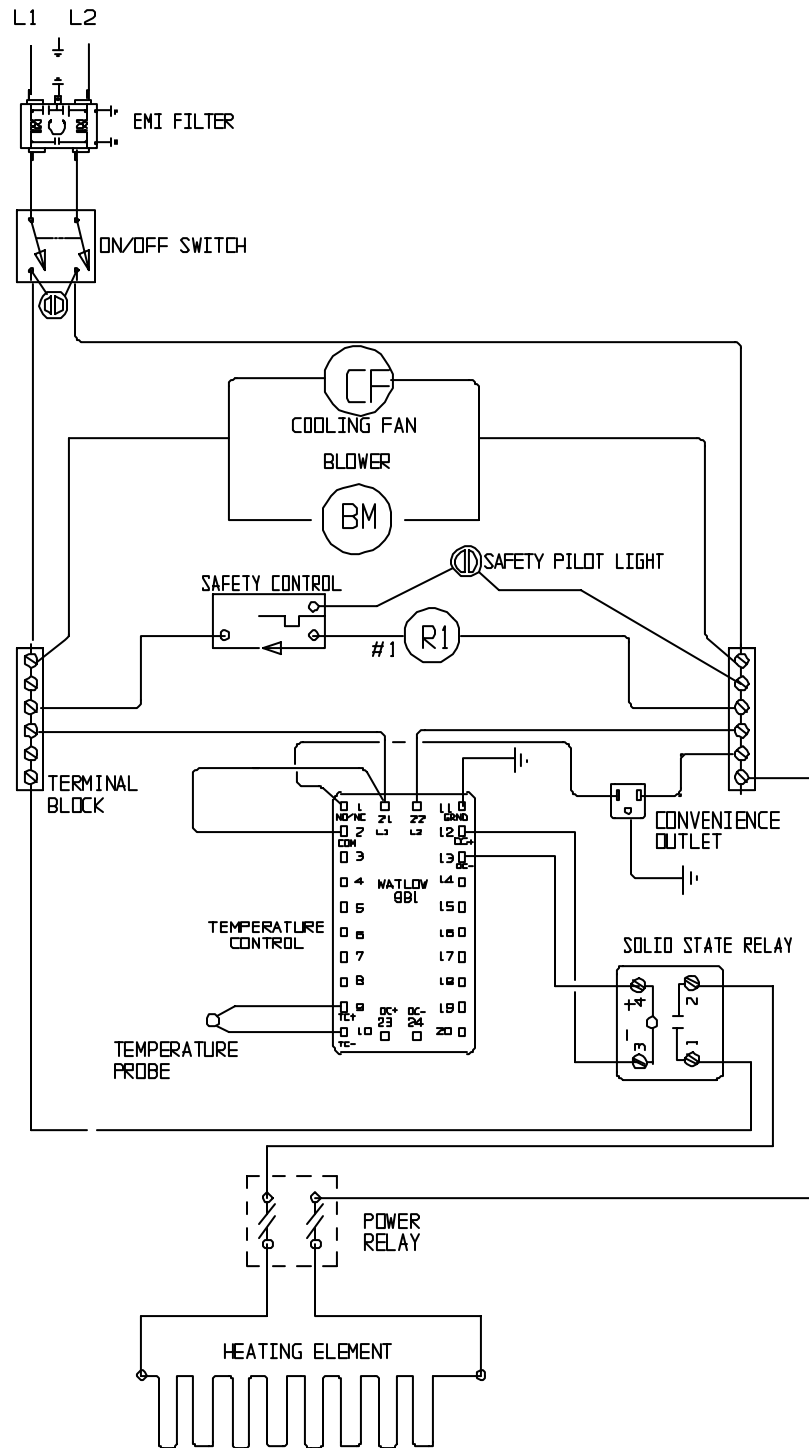

Chapter 6 - Warranty

TestEquity Inc. Limited Warranty

TestEquity Inc. (TestEquity) warrants Environmental Chambers (Equipment) manufactured by TestEquity and supplied under this contract to be free from defects in materials and workmanship under normal use and proper maintenance.

TestEquity will repair or replace any defective part for a period of THREE YEARS from the date of invoice. TestEquity reserves the right to require any defective part be returned, freight prepaid, to TestEquity's factory or to inspect any defective part at the Purchaser's site. TestEquity shall have sole discretion to determine whether any part is defective and whether any defective part will be repaired or replaced. This limited warranty shall extend to any standard chamber accessory and component part which is normally sold by TestEquity. Non-standard accessories and component parts specified by the Purchaser shall be warranted only to the extent of the original manufacturer's warranty, if any exists.

If the repair or replacement is performed in the FIRST YEAR from the date of invoice, TestEquity will also pay for the labor associated with the repair at the Purchaser's site, subject to TestEquity's prior approval. During the SECOND and THIRD YEAR of the warranty period, Purchaser will be responsible for the installation and cost of installation of replacement or repaired parts. Purchaser shall notify TestEquity in writing of any alleged defect within 10 days after its discovery within the warranty period. TestEquity reserves the right to satisfy the labor portion of this limited warranty either through its own service personnel or an authorized agent. In order to provide expeditious service, TestEquity reserves the right to satisfy its limited warranty obligation by sending replacement parts to be installed by the Purchaser if they can be installed easily without special tools or training. TestEquity reserves the right to satisfy this limited warranty by requiring the Purchaser to return the Equipment to TestEquity when such return is feasible.

The following parts are excluded from this limited warranty and are sold as-is or are considered expendable: interior light bulb, viewing window, paint and cosmetic surface finishes and treatments, and port plugs.

This limited warranty shall extend in full to Equipment installed within continental United States and Canada. For all other locations, Purchaser is responsible for all labor costs for repairs or parts installation, and for all shipping costs associated with providing replacement parts.

This limited warranty does not cover: (1) Defects or damages arising as the result of shipment by common carriers or private transportation, unless TestEquity undertakes shipment and transportation of the Equipment to Purchaser's site or contractually assumes the risk of damage to the Equipment in shipment; (2) Defects or damages arising out of, or as the result, of mishandling, modification, or improper start up, installation or maintenance of the Equipment (including start up, installation or maintenance not in accordance with TestEquity's written procedures); (3) Defects or damages resulting from, or arising out of, abuse, misuse, neglect, intentional damage, accident, fire, flood, earthquake, or any other act of God.

This warranty as to Equipment is LIMITED to repair or replacement of parts or Equipment in the determination of TestEquity Inc. THE FORGOING LIMITED WARRANTY IS IN LIEU OF ALL OTHER WARRANTIES INCLUDING THE IMPLIED WARRANTIES OF FITNESS FOR A PARTICULAR PURPOSE AND MERCHANTABILITY. TestEquity Inc. DISCLAIMS ANY LIABILITY FOR ANY DAMAGES RESULTING FROM DELAY OR LOSS OF USE IN SERVICE OR REPAIR, OR FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES ARISING OUT OF OR IN CONNECTION WITH THE USE OR PERFORMANCE OF THE EQUIPMENT, EXCEPT AS STATED IN THIS PARAGRAPH.

This limited warranty cannot be modified in any way except in writing by both TestEquity and Purchaser. Invalidation of any one or more of the provisions of this limited warranty shall in no way affect any of the other provisions hereof, which remain in full force and effect.

This limited warranty shall be extended only to the first Purchaser of this Equipment and is not transferable.